


INSTRUCTION MANUAL
CAT. 64401-10, 64401-50 Tall Slanted Shelf Cart




Electron Microscopy Sciences
1560 Industry Road Hatfield, PA 19440
TEL: 215-412-8400 FAX: 215-412-8450 TOLL FREE: 1-800-523-5874
EMAIL: sgkcck@aol.com WEB: www.emsdiasum.com

Thank you for choosing Electron Microscopy Sciences for your storage and organizational needs! This guide will show you step by step how to assemble your Tall Slanted Shelf Cart.


INCLUDED COMPONENTS

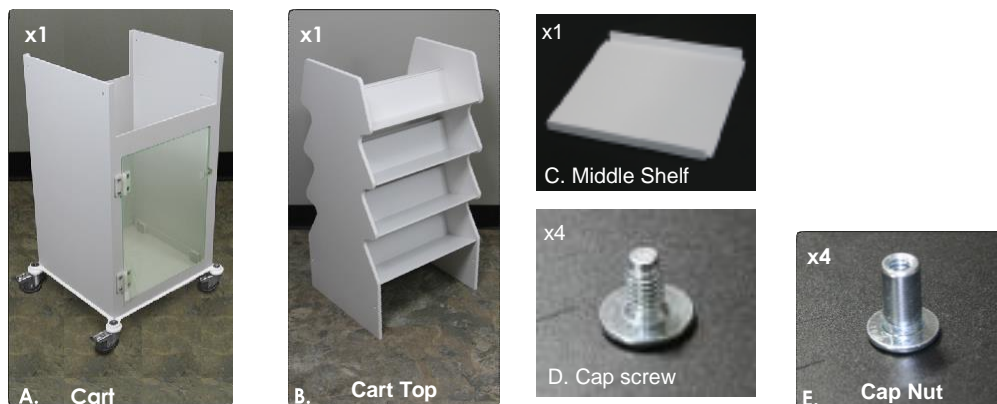

UNPACKING THE CART

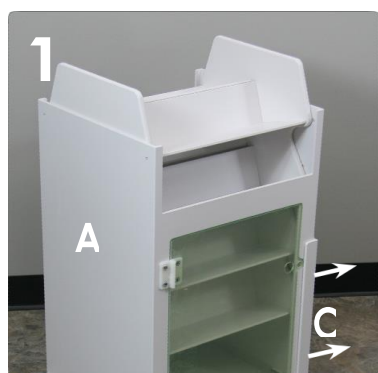


The main pieces of the cart come nested together. Remove the middle shelf "C" from the right side of the cart base "A" and set aside.

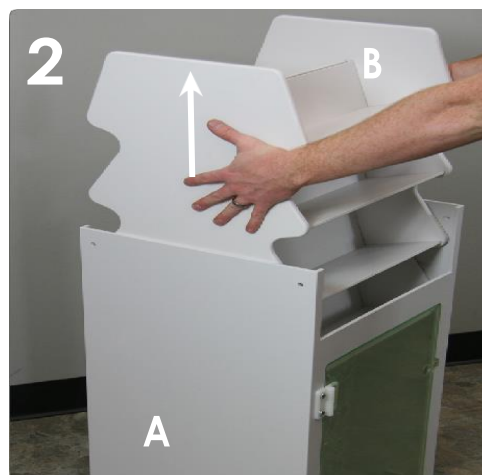


Lift the cart top "B" out from inside the cart base "A".


Electron Microscopy Sciences
 1560 Industry Road Hatfield, PA 19440
 TEL: 215-412-8400 FAX: 215-412-8450 TOLL FREE: 1-800-523-5874
 EMAIL: sgkcck@aol.com WEB: www.emsdiasum.com

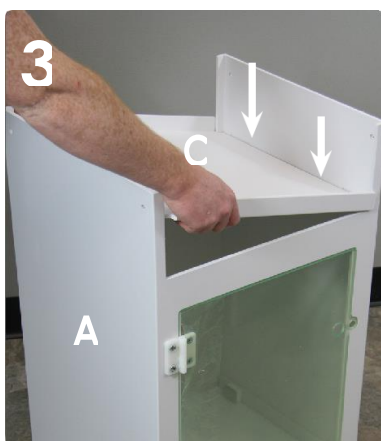

3
Lay the middle shelf "C" with long tabs facing up into the grooves on the inside of the cart base "A". They will pop into place.

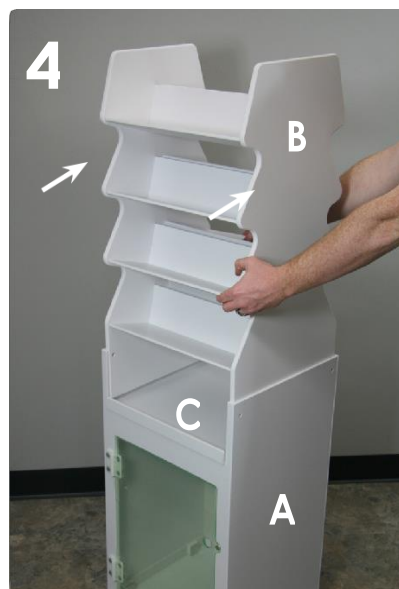

4
Next place the cart top "B" inside the cart base "A". The cart top should rest on top of the middle shelf "C" installed in the last step.

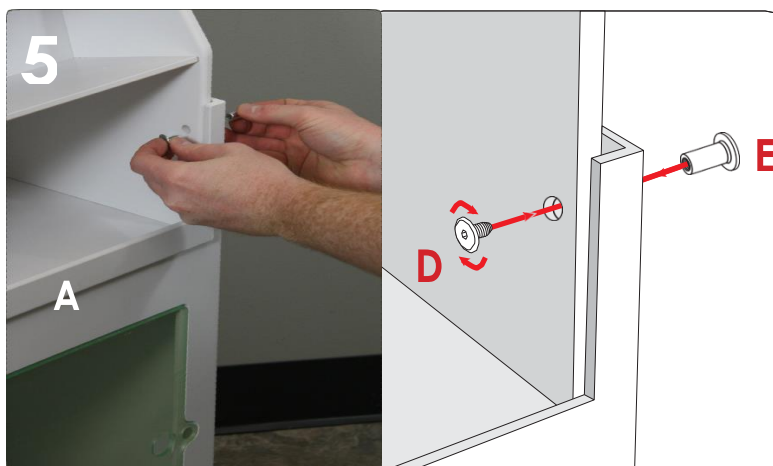

5
Insert the cap nut "E" through the holes in the cart base "A" and cart top "B" as shown above. Tighten the cap screw "D" into the cap nut "E". Repeat with all four holes.

Electron Microscopy Sciences
1560 Industry Road Hatfield, PA 19440
TEL: 215-412-8400 FAX: 215-412-8450 TOLL FREE: 1-800-523-5874
EMAIL: sgkcck@aol.com WEB: www.emsdiasum.com