


INSTRUCTION MANUAL

CAT. 24145-10, 24145-25 Red-Z – Fluid Control Solidifier



Electron Microscopy Sciences
1560 Industry Road Hatfield, PA 19440
TEL: 215-412-8400 FAX: 215-412-8450 TOLL FREE: 1-800-523-5874
EMAIL: sgkcck@aol.com WEB: www.emsdiasum.com

With the Red Z – Fluid Control Solidifier, you can solidify potentially infectious blood and body fluid spills. It is fast-acting and easy to use, eliminates splashing by solidifying and turning spills, deodorizes spills, and is a cost effective cleanup solution.

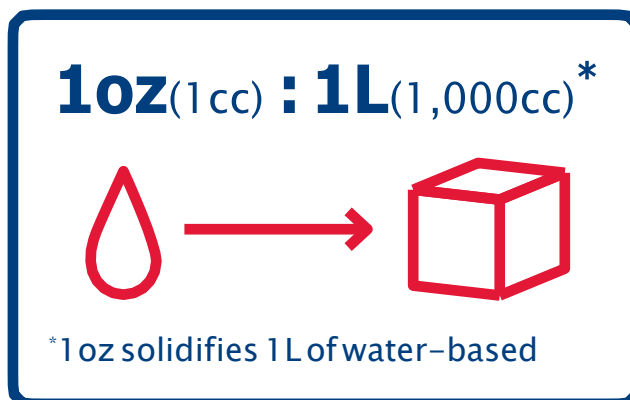

DIRECTIONS

1. Wear appropriate protective clothing.
2. Sprinkle Red-Z Fluid Control Solidifier evenly over spill.
3. Allow sufficient time for solidification to occur.
4. Remove solidified material with scoop and scraper.
5. Clean and disinfect contaminated areas after use and wash thoroughly.



Electron Microscopy Sciences
1560 Industry Road Hatfield, PA 19440
TEL: 215-412-8400 FAX: 215-412-8450 TOLL FREE: 1-800-523-5874
EMAIL: sgkcck@aol.com WEB: www.emsdiasum.com