


INSTRUCTION MANUAL
CAT. 72359-10 Mini Microcentrifuge




Electron Microscopy Sciences
1560 Industry Road Hatfield, PA 19440
TEL: 215-412-8400 FAX: 215-412-8450 TOLL FREE: 1-800-523-5874
EMAIL: sgkcck@aol.com
WEB: www.emsdiasum.com

Operating procedures:


- After unpacking, place the Mini Microcentrifuge on a flat, clear surface
- To open, press firmly down on release tab while lifting dome
- Insert tubes into rotor completely. Be sure that each tube is securely closed and of equal volume. If necessary, add dummy-tubes to balance.
- Load rotor as indicated below. Properly balanced loads are essential for safe and quiet operation.

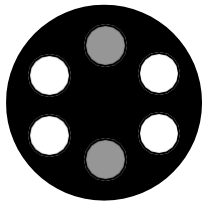


2 tubes

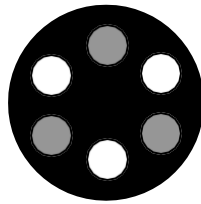

3 tubes

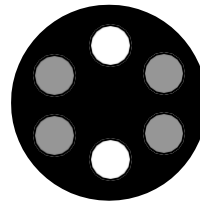

4 tubes

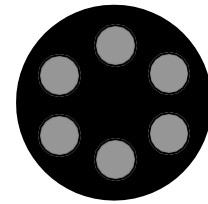

6 tubes

- Close dome securely and activate side switch. Device will not operate with dome open.
- To stop device, deactivate side switch. Wait until device has completely stopped to release dome. Remove samples.

Safeguard

- Do not immerse device in water. Injury due to shock or fire may result.
- Do not operate device with dome open, or open dome before rotor has ceased spinning.
- Do not insert non-standard tubes or other foreign articles into rotor.
- Do not operate under uneven loads. Destabilized operation may result in damage to the device or injury to the user.
- Store item at room temperature in a dry area. Do not expose to sun, moisture or extreme temperature for prolonged periods of time.